

## SpeedVac

- **Preheating:**  
15min before starting the run.  
Set to „medium“ to reach 43°C or to „high“ for 65°C. („low“ means room temperature!)
- Start refrigerated vapor trap (on floor). Let it cool down to -85°C. Then vacuum pump will start automatically when this temperature will be reached.
- Load samples symmetrically, close the cover and set the concentrator switch to „on“.
- Allow rotor to reach full speed (ca. 20sec)
- Continue run until a desired time has elapsed or until predetermined vapor pressure is reached.
- Shut off the speedVac concentrator. The vacuum will be released, the lid can be opened and the samples can be taken out.
- Shut off the refrigerated vapor trap. The vacuum pump will also turn off.